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Please read the Instruction Manual carefully, for correctly grasping the installation and using method of this product. Please keep properly this Instruction Manual after reading, for the usage during troubleshooting and maintenance in the future.

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Tel: 86-0532-66993768

E-mail: [cmtech@sina.com](mailto:cmtech@sina.com)

Address: No. 7 Wenzhou Road, City Northern District, Qingdao City, China

Production base: No. 3 Xinghai Road, Liuting Street, Chengyang District, Qingdao

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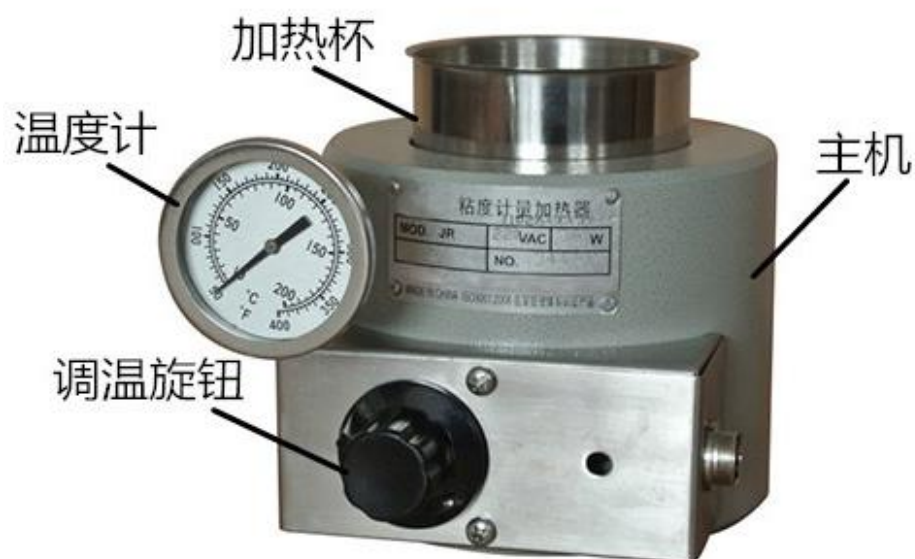
## I. Overview

Viscosity metering heater is used with our company's model:1100,1103,1104 and 6-speed rotating viscometer, it is used for measuring the rheological parameter by simulating the constant temperature of drilling fluid under the well, it also can be used for carrying out the heating and constant temperature for other fluid. The instrument is used widely in petroleum, geological exploration and others.

## II. Model and Specifications

Model: 1106

## III. Main Technical Parameters of Instrument

Name	Technical parameters
Rated power source	220V±5% 50/60Hz
Rated power	300W
Control temperature	Control the room temperature to 93℃
Time of cup warming	Less than 30 minutes

## IV. The Instrument Structure

## **V. Instrument Operation**

1. Connect the power source, adjust the thermostat knob to zero.
2. Insert the bimetal thermometer into the thermometer jack on the left side of heater, rotate the thermostat knob to let the heating rod starts to work.
3. Observe the thermometer reading. When the reading is high than the required temperature around 10°C, adjust the knob to stop the warming, the heater starts to keep the constant temperature.
4. Pour the sample into the sample cup, and then put the sample cup into the heater.
5. Put the bimetal thermometer into the drilling fluid, observe the thermometer reading, when the reading reaches to the required temperature, adjust the knob, the sample cup starts to keep the constant temperature.
6. Place the heater into the positioning hole of the tray of six-speed rotating viscometer, so as to carry out the viscosity test.

## **VI. Instrument Care and Maintenance**

1. Clear and dry all the components and parts for the stand-by usage, place the instrument in the dry environment.
2. When moving, maintaining and making the service for the instrument, please take up and place it gently, so as to avoid the components are damaged and the accuracy and usage are influenced.
3. If the sample liquid or water splashes on the instrument, wipe it gently in time, so as to avoid the equipment is damaged and causes the accident.

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©版权所有 青岛创梦仪器有限公司
Qingdao Chuangmeng Instrument Co., Ltd. Packing list
Manufacturing enterprise: Qingdao Chuangmeng Instrument Co.,Ltd.