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© all rights reserved, Qingdao Chuangmeng Instrument Co., Ltd

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Please read the Instruction Manual carefully, for correctly grasping the installation and using method of this product. Please keep properly this Instruction Manual after reading, for the usage during troubleshooting and maintenance in the future.

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Address: No. 7 Wenzhou Road, City Northern District, Qingdao City, China

Production base: No. 3 Xinghai Road, Liuting Street, Chengyang District, Qingdao

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The mixer produced by Qingdao Chuangmeng Instrument Co., Ltd. It is mainly used for the configuration of drilling fluid base slurry. The mixer has the advantages of simple structure, reasonable design, reliable quality, high starting torque and larger stirring force, convenient speed regulation, free adjustable angle, adjustable position, flexible and convenient use and so on. It can be applied to all kinds of drilling fluid laboratory and field, and can be used in pharmaceutical, electronic and scientific research units to prepare the analytical fluid agitation.

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- 1) Connect the 220V power.
- 2) Press the knob to start. The equipment started to run. Rotate the knob, adjust the speed. The rotation speed is increased to the right. The rotating speed to the left is reduced.
- 3) During operation, press the knob to stop the device running.

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- 1、 Clean all parts and dry. The instrument is placed in a dry environment.
- 2、 When moving, repairing, or maintaining the instrument. Be careful so as not to cause parts distortion, affect accuracy and use.
- 3、 The mixing shaft is not collided. So as not to bend the deformation and damage the instrument.
- 4、 The motor and governor should be checked regularly. It is forbidden to put corrosive chemicals together with mixers.

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Production address: No. 3 Xinghai Road, Liuting Street, Chengyang District, Qingdao

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**Qingdao Chuangmeng Instrument Co., Ltd. Packing list**
Manufacturing enterprise: Qingdao Chuangmeng Instrument Co.,Ltd.